

AN

Sterile Diamine, Disinfectant Cleaner

Low residue Quat free disinfectant / cleaner. Broad spectrum, effective against bacteria, fungi and viruses. TB effective.



InSpec AN for hard surface disinfection

GENERAL GUIDANCE

USE INSTRUCTIONS:

InSpec AN is designed for spraying, wiping and mopping applications. Sprays hold approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate bucket for mopping. 100ml concentrate make up to 5L with water for mopping. Apply to surface to ensure complete coverage.

TECHNICAL DATA:

Appearance: Straw coloured clear liquid free from insolubles, pH10.6 approximately. Data is typical of normal production and should not be taken as a specification.

STORAGE:

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is provided in a separate Safety Data Sheet.

SAFE HANDLING:

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

MATERIAL COMPATIBILITY:

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

BENEFITS

- Low residue
- Good cleaner and disinfectant
- Bactericidal and Yeasticida.
- 1L and 5L formats diluted with WFI quality water
- Defined Shelf life
- Shows virucidal activity Influenza A (H1N1), BVDV, Vaccinia, HBV, HIV
- Defined Shelf life
- Validated In-Use Shelf Life. Contents protected from "suck back" in 1 Litre trigger spray formats
- Multiple bagged
- Quat free ideal rotational partner for InSpec QT and InSpec OX

Microbiological* Minimum Efficacy Testing Contact

EN - 1276 Bacteria 5 mins
EN - 1650 Fungi 15 mins

*Microbiologique / Mikrobiologische / Microbiologica / Microbiológica / Microbiologische

TECHNICAL USER INFORMATION

AN is a sterile disinfectant in pre-diluted and concentrated formats.

Prediluted formats use WFI quality water.

InSpec AN is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped within an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (at 25kGy) to give a sterility assurance level SAL of 10⁻⁶.

InSpec AN comes with key certification such as irradiation, sterility, endotoxin and analysis.

Trigger spray has 4 month in-use shelf life.

For screw cap formats use the entire contents after opening.

InSpec AN PRODUCTS

PRODUCT	PACK SIZE	CARTON SIZE	PRODUCT CODE
InSpec AN 1L Trigger	1 litre trigger spray	6 x 1L	ANWFI20-900MLS
InSpec AN 5L Screw Cap	5L screw cap	2 x 5L	ANWFI20-5LS
InSpec AN Unit dose concentrate	100ml	50 x 100ml	ANCNC20-100